



USER MANUAL
AUTOCLAVE CONTROLLER (PRINTER)
THREE VALVE
MODEL NO: (SAC-203)

OPERATION

1. SET Key : Use this key to set parameters

Pressing the **SET** Key will scroll through following parameters.

- SET JACKET TEMPERATURE Set point for temperature of Jacket.
- SET CHAMBER TEMPERATURE Set point for temperature of chamber.
- STERILIZATION TIME: Time, for which Chamber temp. is keep at set point.
- DRY TIME: Time, for which dry cycle has to be run after soaking.
- PRINT TIME: Time interval for Data Printing.
- SET HIGH TEMPERATURE Set Temperature
- ALARM BAND Temperature Deviation from Set Point to On Alarm.
- CHAMBER PURGING START Start temperature for Purging.
- CHAMBER PURGING STOP Stop temperature for Purging.
- JACKET PURGING START Start temperature for Purging.
- JACKET PURGING STOP Stop temperature for Purging.
- JACKET PURGING TIME Time to open jacket release valve during Purging.
- CYCLE TIME Cyclic On-Off Valve During Purging.
- SECOND CYCLE VALVE ON/OFF Yes or No Selection For Jacket Release after Cycle End
- CHAMBER CONTROL ACTION On-Off
- HYS1 Hysteresis For Chamber Transfer Control Valve.
- JACKET CONTROL ACTION On-Off
- HYS2 Hysteresis For Heater.
- PRESSURE DP: Decimal position for pressure reading.
- PRESSURE RANGE: Range of reading for pressure measurement.
- PRESSURE INPUT: For enabling pressure sensor.
- CHAMBER ERROR
- JACKET ERROR
- SET MINUTE
- SET HOUR
- SET DATE
- SET MONTH
- SET YEAR
- WELCOME MESSAGE

